



MetPrep 1x™
PRECISION GRINDER/POLISHER

*POWERFUL, RELIABLE & DURABLE PRECISION
POLISHER FOR MANUAL SAMPLE PREPARATION*

Quality Products for Metallographic Sample Preparation & Analysis



The **MetPrep 1x™** precision grinding/polishing machine is excellent for manual preparation for standard applications, or when polishing delicate samples using handheld tools with lapping films that require extremely low RPM and consistent low-end torque. The 7" color LCD touchscreen with keypad entry is used to control all functions with an intuitive interface optimized for efficiency. Five (5) user-defined preset buttons allow quick and easy RPM selections from 5-350 RPM in either clockwise or counterclockwise rotation. The timer function allows either count-up or countdown operation.

Durable: Engineered and constructed using precision-machined hard-anodized aluminum and stainless steel components to provide maximum longevity.


Easy to use: Straightforward controls and user-friendly functions for quick, repeatable operation.

Either precision, standard or magnetic 8"/203 mm or 10"/254 mm platens may be used with any plain/adhesive backed or magnetic system disc.

Precision/standard/magnetic platens, accessories and consumables are sold separately.

Item	Description
5-2350	MetPrep 1x™ , 100-240 V, 5-350 RPM Dims: 15" W x 26" D x 9" H (381 x 660 x 230 mm) Weight: 66 lb. (30 kg)

Features:

- Variable platen speed: 5-350 RPM (in increments of 5 RPM) with five (5) adjustable preset buttons
- 7" color LCD touchscreen with keypad entry to control all functions
- Intuitive interface optimized for efficiency and function
- Powerful, quiet 0.7 HP (535 W) Brushless DC motor that provides constant high-torque output throughout the speed range
- Sturdy RIM, aluminum and stainless steel construction
- Clockwise/counterclockwise platen rotation
- Timer for count-up or countdown operation
- Molded one-piece corrosion and impact resistant enclosure
- Overflow diversion to protect internal components
- Quick-change platen design, anodized to resist wear and corrosion
- Low profile, tabletop design
- Electronic coolant control with adjustable flow control valve
- Retractable coolant nozzle that allows quick and easy sample/bowl cleaning
- CE compliant for EU
- Two (2) year warranty
- Designed and manufactured by Allied in the USA 

Retractable Nozzle



Retractable coolant nozzle allows quick and easy sample/bowl cleaning.

LCD Touchscreen



The 7" color LCD touchscreen is used to control all functions and is extremely easy to navigate, allowing greater efficiency among users.

Standard System

Precision/Standard Platens



Precision platens are lapped parallel within 2 microns. Standard aluminum platens are lapped parallel within 25 microns. Both are then hard anodized for maximum wear resistance. They are used with plain/adhesive backed discs, double sided adhesive discs and disc holding bands, or when converting to a magnetic platen using magnetic bases.

Item	Description
5-2005	Standard Platen, 8" (203 mm)
5-6010	Standard Platen, 10" (254 mm)
10-1005	Precision Platen, 8" (203 mm)
10-1007	Precision Platen, 10" (254 mm)


Disc Holding Bands



Disc holding bands are made of anodized aluminum and secure plain back discs to a standard platen. A protrusion around the top of the band allows variable sized discs (± 3 mm) to be secured.

Item	Description
125-10011	8" (203 mm)
125-10012	10" (254 mm)

Double Sided Adhesive Discs (Pk/10)

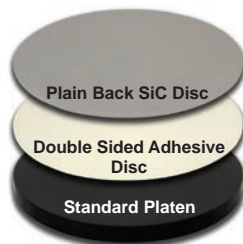


Double sided adhesive discs are adhered to a standard platen and used to secure plain back silicon carbide abrasive discs. Each provides 25-75 disc pulls, depending on grit size, cycle time and force.

Item	Description
52-10005	8" (203 mm)
52-10003	10" (254 mm)

Magnetic System

Magnetic Platens

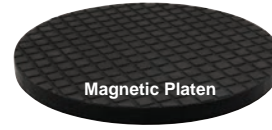

Magnetic platens combine a precision or standard aluminum platen with a cross-hatched magnetic surface for maximum flatness and adherence. Only one platen is needed throughout the grinding and polishing process, allowing all magnetic system cloths/abrasive discs or ferromagnetic discs to be utilized.

Item	Description
5-2005M	Standard Magnetic Platen, 8" (203 mm)
5-6010M	Standard Magnetic Platen, 10" (254 mm)
10-1005M	Precision Magnetic Platen, 8" (203 mm)
10-1007M	Precision Magnetic Platen, 10" (254 mm)


Ferromagnetic Support Discs



These reusable ferromagnetic stainless steel discs feature a nonslip backing and allow adhesive backed cloths/abrasive discs to be changed quickly and easily when using magnetic platens.

Item	Description
55-51000	8" (203 mm)
55-51002	10" (254 mm)

PTFE Ferromagnetic Support Discs

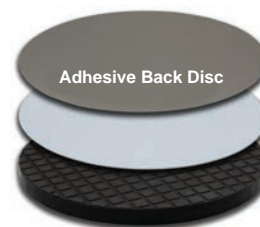

These reusable, durable discs combine a PTFE surface with a ferromagnetic stainless steel substrate. This allows adhesive backed abrasive discs/polishing cloths to be easily removed after use without any adhesive transfer. The nonslip backing keeps the disc in place, even for coarse grinding applications.

Item	Description
55-52000	8" (203 mm)
55-52002	10" (254 mm)

Global Distribution of The World's Finest Products



Brochure #: MP1x-2021

Equipment is subject to change due to innovations/improvements at any time.



Designed and Manufactured by
Allied High Tech Products, Inc.
in the USA

ALLIED
HIGH TECH PRODUCTS, INC.

2376 E. Pacifica Place, Rancho Dominguez, CA 90220
310.635.2466 Worldwide / 800.675.1118 U.S. & Canada
Visit us online! AlliedHighTech.com